

# Software Release Note

## SIMS Processing

Release 2.6 Patch #3

### Release Date

Date	Release	Applicability
December 2, 2024	2.6.3	This software release applies to <ul style="list-style-type: none"><li>• Corrections R1760</li></ul>

## Modifications in Software Release 2.6.3 for SIMS Processing

### 1. List of Bugs Fixed

ID	Application	Description
1760	WinCurve	Fixes a bug relating to the range of 'constant as reference' signal. For quantified profiles, when returning to a c/s profile to reposition the range of the "constant as reference" signal, the reference signal was applied to the first profile displayed in the graph window and not the selected profile. This has been corrected. The range of 'constant as reference signal' is now applied to the selected profile.