

Software Release Note AKONIS Pilot

Release 1.0 Patch #12

Release Date

| Date | Release | Applicability |
|-------------------|---------|---|
| November 28, 2024 | 1.0.12 | This software release applies to <ul style="list-style-type: none">• Corrections : R1760, R1776 |

Modifications in Software Release 1.0.12 for AKONIS Pilot

1. List of Bugs Fixed

| ID | Application | Description |
|------|-------------|---|
| 1760 | WinCurve | Fixes a bug relating to the range of 'constant as reference' signal. For quantified profiles, when returning to a c/s profile to reposition the range of the "constant as reference" signal, the reference signal was applied to the first profile displayed in the graph window and not the selected profile. This has been corrected. The range of 'constant as reference signal' is now applied to the selected profile. |
| 1776 | Vacuum | Fixes a bug relating to the opening of the recorder window. The recorder window now opens correctly. |